



Changsha YiGuang Pharmaceutical Machinery Co.,Ltd

Address: 2/F, Building A1, Lugu International Industrial Area, Yuelu District, Changsha, Hunan, China
Tel:86-731-85634411 Fax:86-731-85634411 Email:felicia@csygyj.com www.labpharmachine.com

SHB Three-dimensional mixer series



Introduction:

SHB series laboratory three-dimensional motion mixer is a novel material mixer widely used in pharmaceutical, chemical, food, light industry and other industries and scientific research units. This

Attn:Felicia
Mobile:+86-15013509826
Skype:felicia@csygyj.com
Wechat:ambostrade

Title:Sales Manager
Email:felicia@csygyj.com
Whatspp/Viber:+86-15013509826



Changsha YiGuang Pharmaceutical Machinery Co.,Ltd

Address: 2/F, Building A1, Lugu International Industrial Area, Yuelu District, Changsha, Hunan, China
Tel:86-731-85634411 Fax:86-731-85634411 Email:felicia@csygyj.com www.labpharmachine.com

machine can mix powder or granular materials with good fluidity very uniformly materials. When the machine is working, due to the multi-directional operation of the mixing barrel, the flow and diffusion of various materials are accelerated during the mixing process. At the same time, it avoids the segregation and accumulation of the specific gravity of the materials caused by the centrifugal force of the general mixer. Good mixing effect can be achieved.

Working Principle:

Driven by the drive shaft, the loaded cylinder makes repeated translational, rotating, and tumbling combined motions, prompting the material to perform circular, radial, and axial three-way compound movements along the cylinder, thereby realizing a variety of materials. Mutual flow, diffusion, accumulation, doping, to achieve the purpose of uniform mixing.

Technical Parameter:

Attn:Felicia
Mobile:+86-15013509826
Skype:felicia@csygyj.com
Wechat:ambostrade

Title:Sales Manager
Email:felicia@csygyj.com
Whatspp/Viber:+86-15013509826



Changsha YiGuang Pharmaceutical Machinery Co.,Ltd

Address: 2/F, Building A1, Lugu International Industrial Area, Yuelu District, Changsha, Hunan, China
Tel:86-731-85634411 Fax:86-731-85634411 Email:felicia@csygyj.com www.labpharmachine.com

Model number	Mixing tank volume (L)	Maximum loading volume (L)	Spindle speed (r/min)	Power (kw)	Dimension (mm) (L×W×H)	Weight (kg)
SHB-20	20	16	0-30 Frequency Control	0.55	740X1000X850	120
SHB-50	50	40	0-30 Frequency Control	1.5	1000X1400X1300	320
SHB-100	100	80	0-12 Frequency Control	2.2	1200X1550X1400	500
SHB-200	200	160	0-12 Frequency Control	4	1300X1600X1500	800
SHB-300	300	240	0-12 Frequency Control	4	1350X1650X1550	1000
SHB-400	400	320	0-12 Frequency Control	4	1500X2200X1500	1200
SHB-600	600	480	0-12 Frequency Control	5.5	1850X2500X1750	1500
SHB-800	800	640	0-12 Frequency Control	7.5	2100X2650X2000	1700
SHB-1000	1000	800	0-12 Frequency Control	7.5	2150X2800X2100	1800

Attn:Felicia
Mobile:+86-15013509826
Skype:felicia@csygyj.com
Wechat:ambostrade

Title:Sales Manager
Email:felicia@csygyj.com
Whatspp/Viber:+86-15013509826