



Changsha YiGuang Pharmaceutical Machinery Co.,Ltd

Address: 2/F, Building A1, Lugu International Industrial Area, Yuelu District, Changsha, Hunan, China
Tel:86-731-85634411 Fax:86-731-85634411 Email:felicia@csygyj.com www.labpharmachine.com

V-type mixer series



Attn:Felicia
Mobile:+86-15013509826
Skype:felicia@csygyj.com
Wechat:ambostrade

Title:Sales Manager
Email:felicia@csygyj.com
Whatspp/Viber:+86-15013509826



Changsha YiGuang Pharmaceutical Machinery Co.,Ltd

Address: 2/F, Building A1, Lugu International Industrial Area, Yuelu District, Changsha, Hunan, China
Tel:86-731-85634411 Fax:86-731-85634411 Email:felicia@csygyj.com www.labpharmachine.com

Introduction:

This machine is suitable for mixing dry powder and granular materials of pharmacy, chemical industry and food. The mixing barrel of this machine has a unique structure, uniform mixing, high efficiency and no accumulation of materials. The whole machine is simple in structure and easy to operate. The outer surface and the material contact part are made of high-quality stainless steel, beautiful in appearance, easy to maintain and clean.

Working Principle:

Using the principle that the lengths of the two cylindrical barrels of the mixer are not equal, and the rows are asymmetrical, when the mixer rotates, the materials decompose to combine, and from combination to decomposition. Due to the different planes of the materials, there is also a lateral force, which pushes the materials to communicate each time one rotation of the lateral force causes about 25% of the material to flow from one barrel to another, so that the material is mixed horizontally and radially, and the decomposition

Attn:Felicia
Mobile:+86-15013509826
Skype:felicia@csygyj.com
Wechat:ambostrade

Title:Sales Manager
Email:felicia@csygyj.com
Whatspp/Viber:+86-15013509826



Changsha YiGuang Pharmaceutical Machinery Co.,Ltd

Address: 2/F, Building A1, Lugu International Industrial Area, Yuelu District, Changsha, Hunan, China
Tel:86-731-85634411 Fax:86-731-85634411 Email:felicia@csygyj.com www.labpharmachine.com

and combination proceed with each other, so that the material achieves a uniform mixing effect.

Technical Parameter:

Model number	Barrel volume (L)	Working volume (L)	Rotating speed of barrel (r/min)	Motor Power (KW)	Dimension (L×W×H)	Weight (kg)
V-14	14	5.6	28	0.55	930×375×920	72
V-20	20	8	28	0.55	1000×420×880	80
V-30	30	12	23	0.75	1120×405×1050	80
V-50	50	20	20	1.1	1300×460×1200	165
V-100	100	40	18	1.5	1560×705×1500	180
V-200	200	80	15	1.5	1965×860×1750	300
V-300	300	120	13	2.2	2220×920×1950	520
V-500	500	200	12	2.2	2660×1250×2400	600
V-1000	1000	400	12	6.6	3220×1500×2840	1000

Attn:Felicia
Mobile:+86-15013509826
Skype:felicia@csygyj.com
Wechat:ambostrade

Title:Sales Manager
Email:felicia@csygyj.com
Whatspp/Viber:+86-15013509826